

Applied Biosystems™ 3130/3130xl Genetic Analyzers:

Starting Up the Instrument

This datasheet describes how to start or restart the system after performing a long-term or short-term shutdown of the 3130/3130xl Genetic Analyzer Instrument. Step-by-step instructions for the start-up procedures can be found in the [Applied Biosystems™ 3130/3130xl Genetic Analyzers 3130 Series Data Collection Software 4 Getting Started User Guide \(Cat. No. 4477796\)](#).

Turning on the Instrument

1. Turn on the computer and log in as the local administrator.
2. Turn on the power button on the instrument and wait for the green light.
3. Launch the Data Collection Software and wait until all 4 services turn green.

If you are restarting the system after performing a long-term shutdown, use the reactivation option in the Install Array wizard.

Launching the Wizards

1. Follow the instructions for “Turning on the Instrument”.
2. Under GA Instruments on the left side, click on the Instrument Name (e.g., 3130). The Wizards menu should now appear at the top of the screen.