

Release Notes

SimpliAmp™ Thermal Cycler Firmware v1.0.5

For SimpliAmp™ Thermal Cycler

Contents

OVERVIEW

FEATURES

DOWNLOAD INSTRUCTIONS

UPGRADE INSTRUCTIONS

LEGAL STATEMENTS

Overview

Firmware v1.0.5 for the SimpliAmp™ Thermal Cycler

This firmware upgrade is for SimpliAmp™ Thermal Cycler only:

A24811 SimpliAmp™ Thermal Cycler

Firmware version 1.0.5 will include all enhancements of previous releases (1.0.3) in addition to the new features listed below. This version will not in any way affect the performance of your SimpliAmp™ cycler except for the specific changes outlined below.

New Features/Changes of version 1.0.5

1. Check connection to Remote Server when Remote Service is enabled at Settings->Instrument Settings->Remote Service screen
2. Enable PCR Essentials mobile application to display Run Complete status
3. Update run monitor screen details to display 3 zones of sample temperature during experiment run
4. No access to New Method/Incubate function under Restricted mode 2 security setting

Downloading the Firmware File

Download the attached file onto a USB memory stick (shown in .update format).

1. **Important: For Internet Explorer users**, right click the above “download” link. Click “Save Target As”. Change the file extension to “.update” and click save
2. Save/Transfer the .update file (not in any folder) on your USB stick.
3. Make sure that no other .update files are present on the memory stick.

Updating your SimpliAmp™ instrument

1. Power on the instrument.
2. Insert the USB memory stick containing the firmware update into one of the USB ports in the SimpliAmp™ instrument.
3. On the Home screen, touch 'Settings'.
4. In the Settings menu, touch 'Maintenance & Services'.
5. Touch 'Software Update'.
6. If you have downloaded the firmware properly, the file will appear on the screen.
7. Select the file on the screen, and press 'Upgrade' at the bottom of the screen.
8. Touch 'Yes' to confirm that you want to upgrade.
IMPORTANT! Ensure that you do not turn off the instrument when the upgrade is in progress.
9. When the upgrade process is complete, the instrument will restart automatically. When it has rebooted, you can remove your USB stick.

LEGAL STATEMENTS

Important licensing information:

These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

For Research Use Only. Not for use in diagnostic procedures.

© Copyright 2014 Life Technologies Corporation. All rights reserved.